

For Research Use Only (RUO)
Not for use in diagnostic procedures

● CHROMagar™ StrepB



For isolation and differentiation
of *Streptococcus agalactiae* (GBS)

CHR  **Magar**
The Chromogenic Media Pioneer

● CHROMagar™ StrepB

For Research Use Only (RUO). Not for use in diagnostic procedures.

For isolation and differentiation of *Streptococcus agalactiae* (GBS)


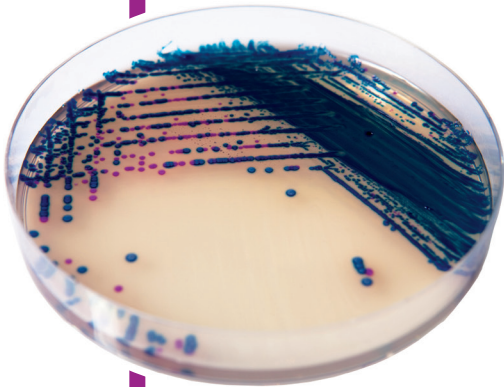


Plate Reading

- Group B *Streptococcus*
→ Mauve
- Other Microorganism
→ blue, colourless or inhibited

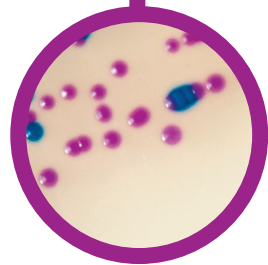

Quality Control Strains

S. agalactiae ATCC® 12386 mauve
S. agalactiae ATCC® 13813 mauve
E. faecalis ATCC® 29212 steel blue
E. coli ATCC® 25922 inhibited
C. albicans ATCC® 10231 inhibited

ATCC® is a registered trademark of the American Type Culture Collection

Medium Description

Powder Base	Total	44.7 g/L
	Agar	15.0
	Peptones and yeast extract	20.0
	Salts	7.5
	Chromogenic mix	2.2
Storage at 15/30°C - pH: 7.3 +/- 0.2		
Shelf Life		2 years
+		
2 Supplements (included in the pack)	1 st : Liquid form.....8ml/L	2 nd : Powder..... 0.25g/L
	Storage at 15/30°C	Storage at 2/8°C
	Shelf Life 2 years	Shelf Life 2 years

Please read carefully the instructions for use (IFU document) available on www.drg-international.com

For US Customers: for Research Use Only

Order References

Please use these product references when contacting DRG:

5000 ml pack SB282 25 L pack SB283-25
 Bulk on request

Manufacturer: CHROMagar
 4 place du 18 juin 1940 75006 Paris - France

Distributed in the US by DRG International
www.DRG-International.com